



Weekly Eyewash Inspection Log

Instructions:

Activate emergency eyewash to ensure proper operation. Water should flow within 1 second of valve release and remain flowing.

Ensure there is sufficient water flow. The stream from both eye pieces should be approximately 6 inches in length and meet in the center of the eyewash nozzle.

Ensure the water is clear. If there are any issues, tag the unit as out of service and notify your safety liaison. A work order to UK PPD shall be submitted for correction as a health and safety issue.

When your annual log sheet is complete, file in your blue Laboratory Safety Manual. Records shall be kept for 3 years.

Date	Initials	Date	Initials	Date	Initials