



## NO BIOHAZARDS

**DO NOT use this autoclave for the decontamination of biological and potentially infectious waste. Sterilization of glassware and liquids may be done in this particular autoclave.**

- ◆ **Never put sealed containers, solvents, volatile or corrosive chemicals, or radioactive materials in the autoclave.**
- ◆ **If the cycle did not meet the appropriate operating conditions or aborted, repeat the autoclave cycle. If the cycle fails again, report the malfunction to the responsible person listed below.**
- ◆ **Before unloading the Autoclave, wait for the pressure gauge to drop to zero with zero time remaining before opening the door.**
- ◆ **Using proper heat resistant PPE, carefully remove autoclaved materials after the cycle is completed and the appropriate waiting time for liquids has been observed.**

### Responsible Person for this Autoclave

Name: \_\_\_\_\_

Phone: \_\_\_\_\_



Research Safety

<https://researchsafety.uky.edu/biological-safety>  
[biosafety@uky.edu](mailto:biosafety@uky.edu)